

MDI single use **Harvester** Funnels do away with the necessity of centrifugation for cell harvests in the lab.

These are designed to clarify up to 2 liter of mammalian cell culture harvests within few minutes as compared to hours by centrifugation.

Construction

MDI **Harvester** Funnels are easy to use vacuum filtration devices which house a large area pleated filter. The filter is made of multilayered progressively finer depth filter media. The special design provides high loading capacity for whole cells and high retention efficiencies for cell fragments, organelles, colloids and lipids to give effective protection to downstream 0.2µm sterilizing filters.

Unique Advantage

- Cell harvest clarification of upto 2 liter within minutes
- Low hold-up volume

Specifications

Funnel Volume: 450 mL and 1 liter

Receiver Bottle Capacity: 500 mL, 1 liter and 2 liter

Funnel Adaptor: 45 mm neck for 500mL and 1 liter bottle
53 mm neck for 2 liter bottle

Effective Filtration Area:

400 cm² for up to 500 mL cell harvest
1200 cm² for up to 2 Liter cell harvest

Materials of Construction

Funnel: Polypropylene

Filter Media: Multilayered Microglassfiber

Filter Components: Polypropylene



For 500 mL

For 1 Liter

For 2 Liter

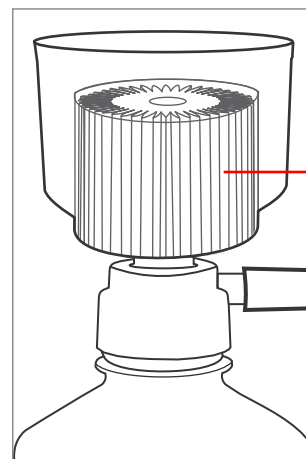


Fig. 1 Large Area Pleated Construction


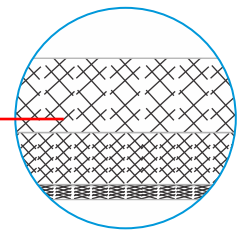


Fig.2 Multilayered Progressive Depth Filtration

Filtration of HEK and CHO Cell Culture

Harvester Funnels were used at customer site for clarification of Human Embryo Kidney (HEK) and Chinese Hamster Ovary (CHO) cell culture. The filtrates were subsequently filtered through 0.2 µm sterilizing grade *AseptiVac KS* vacuum filtration units. The data is shown below:

Cell	Cell Density	Cell Viability	Filtered Volume
HEK	5.2 X 10 ⁶ cells/ml	90 %	1000 mL
CHO	6 X 10 ⁶ cells/ml	80 %	700 mL

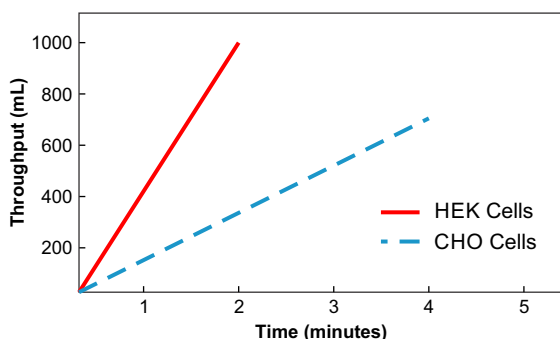


Primary clarification: MDI 500 mL Harvester Funnels (with 1 Liter receiver bottle)

HEK cell culture filtered through the **Harvester** Funnel in 2 minutes and it took 4 minutes to filter the CHO cell culture. The filtrate from HEK cell culture was quite clear. However some cell debris was observed in case of CHO cell culture after filtering 600mL.

Ordering Information

Non Sterile *Harvester* Funnels

- with 500 ml receiver bottle : FHXG5238X05X108
- with 1 liter receiver bottle : FHXG5338X10X108
- with 2 liter receiver bottle : FHXG5338X20X108

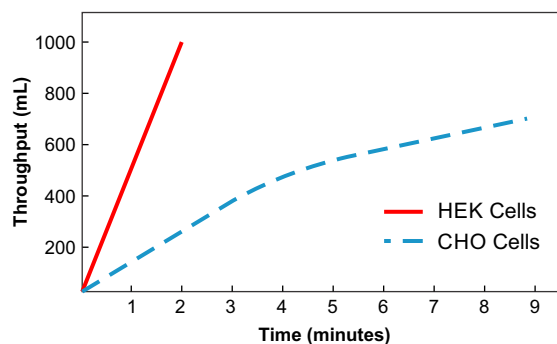

Sterile Filtration (*AseptiVac KS*)

Sterile filtration of the HEK cell harvest filtrate (post primary clarification) through 0.2µm *AseptiVac KS* took only 2 minutes for 1000 mL and took 9 minutes in case of CHO cells for 700mL.

The difference in filtration time was due to the contamination profile of the two cell culture. HEK cell culture with higher cell viability (90%) had much lower cell debris in comparison to CHO cell culture with 80% cell viability.

Important: To minimize frothing, please use the **Harvester Funnel** initially at a low vacuum of <50 mmHg, and increase gradually as required.

Gamma Sterile *Harvester* Funnels

- with 500 ml receiver bottle : FHXG5238X05X308
- with 1 liter receiver bottle : FHXG5338X10X308
- with 2 liter receiver bottle : FHXG5338X20X308

mdi Membrane Technologies INC.

75 Utley Drive STE 103, Camp Hill, PA 17011
 E-mail: rs@mdimembranetech.com
 Phone: +1-717-412-0943, Fax: +1-717-695-9637
 Website: www.mdimembranetech.com

ADVANCED MICRODEVICES PVT. LTD.

20-21, Industrial Area, Ambala Cantt- 133 006, India
 Tel: +91 - 171-2699290, 2699471
 E-mail: info@mdimembrane.com
 Website: www.mdimembrane.com